

LABSTAC



OPERATION MANUAL
STEAM STERILIZER
AUT035-100

www.labstac.com

Index

| | |
|---|----|
| 1. FEATURES..... | 3 |
| 2. TECHNICAL PARAMETERS..... | 4 |
| 3. INSTALLATION REQUIREMENTS..... | 4 |
| 4. Usage Instructions..... | 5 |
| 5. Caution and Maintenance..... | 11 |
| 6. Common Fault Analysis and Troubleshooting..... | 14 |
| 7. PACKING LIST..... | 15 |

BRIEF

The steam sterilizer is the sterilization equipment that can sterilize medical equipment, surgical dressing, medicine liquid, glassware, etc by saturated steam. They enjoy the following merits such as fashionable and attractive appearance, reasonable frame, multiple functions, fast heating, and complete disinfection.

Purpose: they are used to sterilize following goods by medical treatment and public health, scientific research organization: medical equipment, labware, culture medium and non-sealed liquid or preparations.

01. FEATURES

1. This is an automatic control steam sterilizer. This is made by shell, barrel, basket with netting on top, and control system.
2. Shapes are square vertical. The sterilization barrel's shell and themselves are made of stainless steel. Sterilization barrel (pressure vessel) and safety accessories conform to GB 150、《special equipment's safety regulation》 and nationality legal pressure vessel provisions by laws and regulations. They must have nameplates and quality certificates that were registered on the legal pressure vessel safety supervision organization.
3. The safety valve is fixed on the vessel's cover. It can release over pressure to ensure equipment's safety when inner pressure exceeds 0.24MPa.
4. During the heating process, the deflation valve, which is controlled by a temperature controller, will discharge cold air automatically. When the temperature on temperature controller exceeds 102°C, the deflation valve will auto-close after several minutes. The equipment goes into auto intermittent deflated function and raises the vessel's steam saturate degree, which can enhance the sterilization effect.
5. Operating front panel includes a pressure gauge that shows the pressure state clearly.
6. The machine has multiple safety protection. The circuit can automatically shut off when short water in the vessel, heater auto stops operating, protecting the equipment in effect. At the same time, the water indicator light turns red which reminds users to add water until enough.

| Seq. | Volume | Power | Sterilization Time Adjustment Range | Max. Working Pressure | Max. Working Temperature | Net Weight |
|------|--------|-------|-------------------------------------|-----------------------|--------------------------|------------|
| 1 | 35L | 3.5kW | 4-120min | 0.22Mpa | 134°C | 58 kg |
| 2 | 50L | 3.5kW | | | | 80 kg |
| 3 | 75L | 4.5kW | | | | 90 kg |
| 4 | 100L | 4.5kW | | | | 106 kg |

02. TECHNICAL PARAMETERS

1. Power Supply : AC220V \pm 22V , 50Hz \pm 1Hz.
2. Working pressure: 0.22MPa; Working temperature: 116°C to 134°C.
3. During normal operation within the sterilization cycle, the noise should not exceed 65dB (A-weighted).
4. Set sterilization time range: 4 to 120 minutes; Set drying time range: 0 to 240 minutes.
5. Equipment safety class: Class I.
6. Operating environment temperature: +5°C to +40°C; Relative humidity: not more than 85%; Atmospheric pressure 70kPa to 106kPa.
7. The water source used in the sterilizer should not affect the sterilization process or damage the sterilizer and its contents.

03. INSTALLATION REQUIREMENTS

1. Please read the manual carefully before use.
2. According to protection against electric shock, it belongs to class I. No flammable and explosive, anesthetics air or articles exist in installation place.
3. Connect the power source always during the procedure. The sterilizer must keep a 60cm distance from the wall and the height must exceed 1.2m, make sure to operate the equipment safely in (must conform to GB 14048.1 and GB 14048.3 standard) sterilizer's power source. It must have obvious disconnecting device ON-OFF mark (such as special for sterilizer) . The power line must be fixed when it is connected to disconnecting device to prevent from falling off. The power line's diameter is 2.5mm² \times 3. L is the power line. N is the zero line. The yellow-green G is protective grounding. When the procedure is over, please press OFF to close and disconnecting device.
4. Place the equipment on the level and fast floor. Users can connect the power supply correctly according to its nameplate and manual. The equipment must be sure of safety and ground protection (ground resistance shouldn't be more than 2 Ω)
5. Connect the drainage pipe to the drainage port, ensuring that the drainage end of the pipe is not submerged in water to prevent water from being drawn back into the sterilization chamber bucket.
6. For a machine with automatic water inlet function, prepare one bucket (5 gallons) of distilled water (tap water is not recommended, and the water quality should meet the requirements specified in GB 8599 Appendix C). Connect one end of the water inlet pipe to the device's water inlet, and insert the other end into the bottom of the water bucket.

04. Usage Instructions

All documents of the equipment must be supplied to the local inspection organization to register on record.

Sterilizer operators must accept professional training and be familiar with pressure vessel operation according to manual strictly.

Operator must look after the equipment and record the operation state during the sterilizing process, to prevent accidents.

Sterilizers must be checked one time per year at least by an inspection organization. The machine can be used after proving to be qualified.

NOTES :





Please don't put inflammable goods into the sterilizer's barrel.

Please don't put following goods into the barrel: explosive goods, inflammable goods, oxidizing goods, incensive goods, or flammable goods. Otherwise, the equipment will explode, fire or breakdown.


Please don't put the following goods into the barrel such as culture medium that can corrode stainless steel. Otherwise, the equipment will explode, fire, or breakdown.

- Please don't touch the switch or circuit breaker with wet hands, to prevent electric shock or accident.
- Please don't touch the vessel's cover in the sterilizing state as it is high temperature.
- "Please don't open the cover before releasing steam completely": after sterilization, please don't open the cover before the pressure gauge's pointer doesn't return to 0 place. Otherwise, high-temperature steam can burn the operator.

(I) SYMBOLS AND THEIR MEANING


| Symbol | Meaning | Symbol | Meaning |
|---|---------------------|---|---|
| ~ | Alternating current |  | notice ! check and review equipment's documents carefully |
|  | Caution hot |  | Protective grounding |
|  | Test interface | | |

(II) OPERATING STEPS

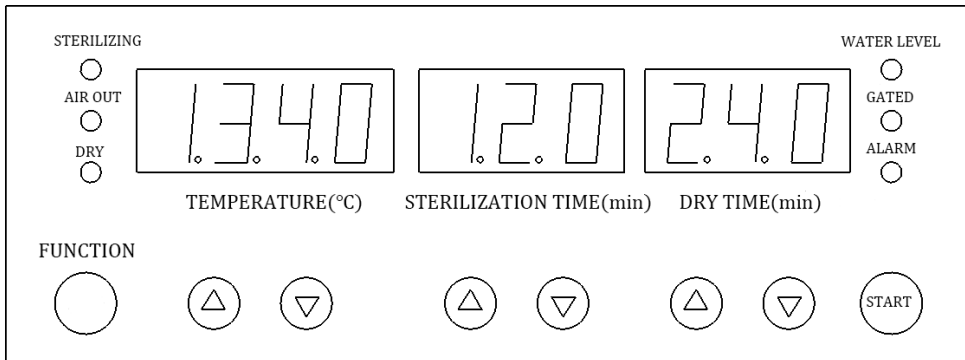



- (1) **Function:** press this button for 5s can choose a function, press one time to display the A model (sterilization function); press again to B mode (sterilization, air release function); press again to display the A model, cycle like that.
- (2) **△ button:** press this button to adjust the temperature display high-temperature setting (range: 50.0~134.0°C), the sterilization time (range: 4 ~ 120 minutes) can increase the set time range. The set time within the drying period (set range: 0 to 240 minutes) can be increased.
- (3) **▽ button:** press this button to adjust the temperature display low-temperature setting (range: 50.0~134.0°C), the sterilization time (range: 4 ~ 120 minutes) can reduce the set time range. The set time within the drying period (set range: 0 to 240 minutes) can be decreased.
- (4) **START:** press this button can open the working function.
- (5) **Water level indicator:** turn red when water is too little in the sterilizer. When the water reaches to a low water level, it turns flashing green. When the water is enough, it displays green light.
- (6) **Gated indicator:** turns off when the sterilizer lid is opened.
- (7) **Alarm indicator light:** turns on when there is an over-temperature alarm or error occurs during work.
- (8) **Temperature display window:** set the state display of set temperature, the working state display current measuring temperature sterilizer.
- (9) **Sterilization display window:** set state display of sterilizing time, display the current surplus of sterilization time working state.
- (10) **Drying Time Display Window:** Displays the set drying time during the setup mode and the remaining drying time during operation.




◆ METHOD OF FUNCTION SETTING

After connecting the power and switching on the circuit breaker, the control panel displays the current temperature in the container and the last run program. Press the "Function" button to select the program operation mode. Then you can use the corresponding  key to set temperature, sterilization time, and drying time respectively. After pressing the "start" button, sterilization indication light will flash (expressed in heating). When the temperature reaches the set value, the sterilization indicator will stay on steadily, and the countdown will begin (indicating the sterilization process). After sterilization, the system will automatically release pressure, and the pressure release indicator will light up. Once pressure release is complete, the drying indicator will light up, and the drying time display window will start counting down (indicating the drying process). The display window will show "END," and a prompt sound will signal the end of the program, allowing the removal of items.

a. Setting Sterilization function

- (1) Press "FUNCTION" button. Display the currently selected working procedure of PROGRAM A, sterilization temperature and sterilization time.
- (2) Set the temperature by pressing the button  which is under the temperature display window.
- (3) Set the sterilization time by pressing the button  which is under the sterilization time display window.
- (4) Press START button after setting well.

b. Setting sterilization, air release, dry function:

- (1) Press the "Function" button. Display the currently selected working procedure PROGRAM B, temperature of sterilization and sterilization time.
- (2) Set the temperature by press the button  which is under the temperature display window.
- (3) Set the sterilization time by press the button  which is under the sterilization time display window.
- (4) Set the drying time by press the button  which is under the drying time display window. (when the drying time is set to 0, there will be no drying after sterilization).
- (4) Press "START" button to start work after setting well.

A. Adding water and Stacking

- (1) Rotate the handwheel counterclockwise for several turns (be cautious to avoid pinching hands and causing injuries) until it reaches the top, lifting the sterilizer cover sufficiently. Push the cover to the right to move the crossbeam and open the sterilizer cover. Remove the sterilization basket.
- (2) Close the water drain valve and pour approximately 5 liters of clean water into the

sterilization drum. The water level should reach the foot of the sterilization drum (below the shelf). For consecutive use, it is essential to replenish water after each sterilization cycle. If there is an automatic water addition function, after function is set, press the start button for automatic water addition. The automatic water addition will continue until the water level indicator light changes from red to a steady green, at which point it will automatically stop.

(3) Properly wrap and arrange the items to be disinfected in the sterilization basket in an orderly manner, leaving gaps between them. This facilitates steam penetration, improving the sterilization effect. When stacking sterilization packages, pay attention to leaving space for the safety valve vent. Otherwise, if the safety valve vent is blocked and cannot release pressure, it may result in the rupture of the sterilizer body.

B. Power On

Connect the power supply that corresponds to the sterilizer label. Set the switch on the control panel to the "On" position. The control panel should display a complete illumination, and a self-check status will be shown for 3 seconds.

C. Sealing

Push the crossbeam towards the left column. Rotate the handwheel clockwise to press the sterilizer cover tightly against the lower flange, exerting force to ensure a complete seal. (When the door control indicator light on the control panel is illuminated, continue rotating until fully sealed. At this point, the following operations can be carried out.)

D. Function Setting

Users should select the appropriate buttons to set the required sterilization temperature and time based on the type of sterilized items.

<Refer to the operation panel instructions. Different items have specific sterilization temperature and time requirements according to the Ministry of Health's "Disinfection Technical Specifications.">

E. Heating

After pressing the start button, the equipment begins operation. The heating indicator on the control panel blinks, indicating the start of heating. The pressure relief valve automatically closes after the temperature reaches 102°C.

F. Sterilization

Once the temperature reaches the set level, the sterilization indicator on the control panel stops blinking (remains solid), indicating the start of sterilization.

G. Pressure Relief (Strictly prohibited when sterilizing glassware and solution culture media)

After sterilization, automatic pressure relief begins (applies to programs in setting mode b). After sterilization, no pressure relief is needed (applies to programs in setting mode a). Allow for natural cooling and pressure reduction.

H. Manual Water Drainage

When necessary, manually open the side drain valve for water drainage. Caution: Be careful to avoid burns during water drainage.

I. Drying

After the water drainage is complete, the automatic drying process will commence. (Note: Drying program is strictly prohibited for liquid sterilization.)

J. Completion

The display window on the panel shows "END," accompanied by a prompt sound, indicating the end. Wait until it cools until the pressure gauge pointer returns to zero, then open the lid and remove the items (be sure to wear protective gloves during retrieval. Due to the temperature drop, extra caution is advised during retrieval. Direct contact with hands may cause burns.) One sterilization cycle is complete. For sterilized glassware and items like solution culture media, open the lid slightly after the pressure gauge returns to "0," wait for 10-15 minutes, and then fully open the lid to remove the sterilized items.

Special Notice: After the work process is completed and "END" is displayed, and when the pressure indicator is zero, you can open the cover to retrieve items. If you find the items too hot and do not wish to retrieve them immediately, it is necessary to manually lift the safety valve to release internal pressure. The cover can be opened later.

Please note: When retrieving items, handle them with care to prevent hot objects from falling and causing burns when opening the door.

During the program execution, if an alarm sounds, the alarm light is on, and the digital display shows a warning symbol (refer to Section VII, Common Fault Analysis and Troubleshooting in this manual for resolution methods), it indicates an error. The errors are as follows:

ERR—1: indicates that the door is not closed.

ERR—2: indicates that the water level is below the low water level.

ERR—3: indicates that the sterilization temperature is too high.

ERR—5: indicates a protection mode for the heating element.

ERR—6: indicates excessive pressure.

ERR—7: indicates an unexpected interruption during the sterilization process.

ERR—8: indicates a manual interruption during the sterilization process.

ERR—9: indicates indoor temperature sensor malfunction.

ERR—10: indicates wall temperature sensor malfunction.

ERR—12: indicates heating film malfunction(for drying function).

ERR—20: indicates a timeout during the heating process.

ERR—21: indicates sterilization failure.

ERR—30: indicates Water inlet failure.

If incorrect settings are noticed after running, press and hold any adjustment key (except "Function" and "Start") for 5 seconds to interrupt the program. Alternatively, turn the switch to the "Off" position, then back to the "On" position, and reselect the desired sterilization temperature and time settings.

NOTE:

For the sterilization temperature correction/sterilization intermittent discharge time and cycle adjustment during the sterilization process: The parameters of this step are related to the equipment working performance. The user should not change the setting easily. If the user does need to adjust the parameters, please adjust as follows :

Temperature Correction Adjustment: Press and hold the start button while turning on the switch. Press the start button once, and the temperature window will flash. Use the plus or minus keys below the temperature window to make corrections. Press the start button again, and the sterilization time window will flash. Use the plus or minus keys below the sterilization time window to adjust the upper limit of the emission gas temperature (105-120°C). Press the start button three or more times, turn off the switch, then turn it on again to complete the correction or adjustment.

Emission Gas Frequency / Intermittent Emission Time and Cycle Adjustment:

For Program A: Hold the function button and the temperature window plus or minus key. Press the start button once, and the temperature window digits will flash. Use the plus or minus keys below the flashing digits to adjust the emission frequency (1-3 times). Press the start button again to adjust the emission time (0-5s). Press the start button once more to adjust the emission cycle (1-60s). Press the start button again to adjust the emission time during gas release (0-10s). Press the start button once more to adjust the emission cycle during gas release (0-10s). Press the start button again to adjust the emission end temperature (110-120°C).

For Program B: Hold the function button and the temperature window plus or minus key. Press the start button once, and the temperature window digits will flash. Use the plus or minus keys below the flashing digits to adjust the emission frequency (1-3 times). Press the start button again to adjust the emission time (0-5s). Press the start button once more to adjust the emission cycle (1-60s).

05. Caution and Maintenance

1. When stacking sterilized items (with the upper cover completely open), the mesh basket should be placed at the center of the barrel (the mesh baskets should be concentric and stacked up and down). It is strictly prohibited to block the vent holes of the safety valve. Sufficient space must be left to ensure smooth ventilation of the equipment. Otherwise, if the vent holes of the safety valve are blocked, it cannot function, leading to accidents.
2. Before each use, check whether the water level in the sterilization barrel is maintained at the specified level.
3. When the sterilizer is in continuous operation and a new sterilization operation is to be performed, allow a 5-minute interval and open the upper cover to give the equipment time to cool down.
4. When sterilizing liquids (only for program A), the liquid container should be placed in a heat-resistant glass bottle, not exceeding 3/4 of its volume. Use a cotton plug for the bottle mouth, and avoid using rubber or cork plugs without holes.
Special note: Do not release steam immediately after the sterilization of liquids. The residual air can only be discharged after the pressure gauge pointer returns to zero.
5. When sterilizing containers such as beakers, flasks, and test tubes, the opening must face downward or be placed horizontally. Placing the opening upward not only makes it difficult to vent but also hinders steam from penetrating the interior, resulting in poor sterilization effectiveness.
6. For items with different types and sterilization requirements, such as dressings and liquids, do not sterilize them together to avoid damage.
7. Confirm sterilization performance: Sterilization performance varies depending on the type of items, quantity, loading method, and container type. Therefore, use sterilization indicators such as an OK card to confirm. Failure to do so may result in accidents or poor sterilization effectiveness.
8. Do not leave sterilized items in the product: This product does not have a storage function for sterilized items. To prevent recontamination, after sterilizing items such as culture media and instruments, promptly remove them and store them in a dedicated safe box.
9. At the end of sterilization (or in case of malfunctions), the pressure gauge pointer must return to zero. Open the manual vent valve (or place the vent valve lever in the vent position) to allow outside air into the sterilizer. Once the vacuum is eliminated, the container lid can be opened (do not approach the lid with your face). When fully opening the lid and removing the mesh basket, be cautious of burns and pinching fingers. (Be sure to wear protective gloves when removing items. Due to the time required for the liquid temperature to decrease, extra care is needed during removal. Direct contact with hands after the operation may cause burns.)
10. After prolonged use, if the pressure gauge shows inaccuracies or cannot return to zero, it should be promptly inspected. Regular checks are recommended (ideally every 6 months) by comparing it with a standard pressure gauge. If abnormalities are found, a new gauge should be replaced. Every six months, apply an appropriate amount of high-temperature grease to the horizontal beam screw.
11. Keep the equipment clean and dry at all times. The rubber sealing gasket may age over time and should be replaced regularly (recommended every 12 months).
12. During operation, if the pressure gauge indicates over 0.24 MPa and the safety valve does not activate, immediately cut off the power, open the manual vent valve, wait for the

pressure gauge to indicate 0 MPa for 1-2 minutes, then open the container lid and promptly replace the safety valve.

13. When troubleshooting requires the replacement of components, it must be done by qualified personnel with professional training or authorized personnel from the manufacturer. Disconnect the power circuit breaker, release the residual air in the container, zero the pressure gauge pointer, and send the safety valve for inspection at least once a year. Lift and lower the handle on the safety valve core several times each week to prevent valve core scaling and air leakage.

14. Do not modify for use: Unauthorized disassembly or repair is strictly prohibited, except for maintenance technicians. Otherwise, injuries may occur due to fire or abnormal operation. If screws or nuts are found to be loose during routine use, tighten them promptly to ensure normal operation.

15. Cleaning (recommended monthly): Empty the water from the container and disconnect the power circuit breaker. After the container has cooled, remove the shelves and mesh basket from the container. Use a brush to remove scale deposits on the inner walls, heating elements, and water level probes. Then, use a towel dipped in alcohol or distilled water to wipe the inner surfaces. After drying, place the shelves and mesh basket back into the container. Store the equipment in a well-ventilated room with non-corrosive gases.

16. The interlock device of the door is composed of a self-locking mechanism and a door control switch. During regular use, ensure the flexible extension and retraction of the self-locking mechanism (recommended quarterly inspection). Do not forcefully open the container lid before the pressure inside the container has been released to zero, as it may damage the self-locking mechanism. The door control switch should maintain flexible contact points. If contact points are stuck or indicator lights are not working properly, prompt maintenance is required.

17. Do not allow children to use or play with the equipment: Otherwise, it may result in burns, electric shock, or injuries.

18. Do not open the drainage valve while the equipment is in operation: When draining the heated water, wait for at least 2 hours after sterilization is complete and ensure that the lid is open before starting the drainage operation. Failure to do so may result in burns or accidents due to the release of high-temperature hot water.

19. Fuse replacement: If the power indicator light does not illuminate after connecting the sterilizer power cord to the power source, disconnect the circuit breaker. Using a Phillips screwdriver, open the back cover of the sterilizer to inspect the fuse (specifications: R015, 10×38, 32A). If the fuse is damaged, replace it immediately.

20. Before moving, transporting, or storing, drain the water from the container, seal the sterilizer lid, disconnect the circuit breaker, and remove the power cord. When moving (release the caster locks first), push the sterilizer on a flat surface. During transportation and storage (secure the sterilizer in the packaging box), prevent loosening and abrasion. Avoid tipping to prevent injuries or accidents.

21. Transportation and storage environmental conditions:


Ambient temperature: -40°C to +55°C

Relative humidity range: 10% to 80%

Atmospheric pressure range: 500 hPa to 1060 hPa

22. At the end of its service life (8 years from the production date indicated on the product label), this product and its accessories should be voluntarily handed over to qualified entities or individuals for recycling and disposal.

23. The wiring diagram:

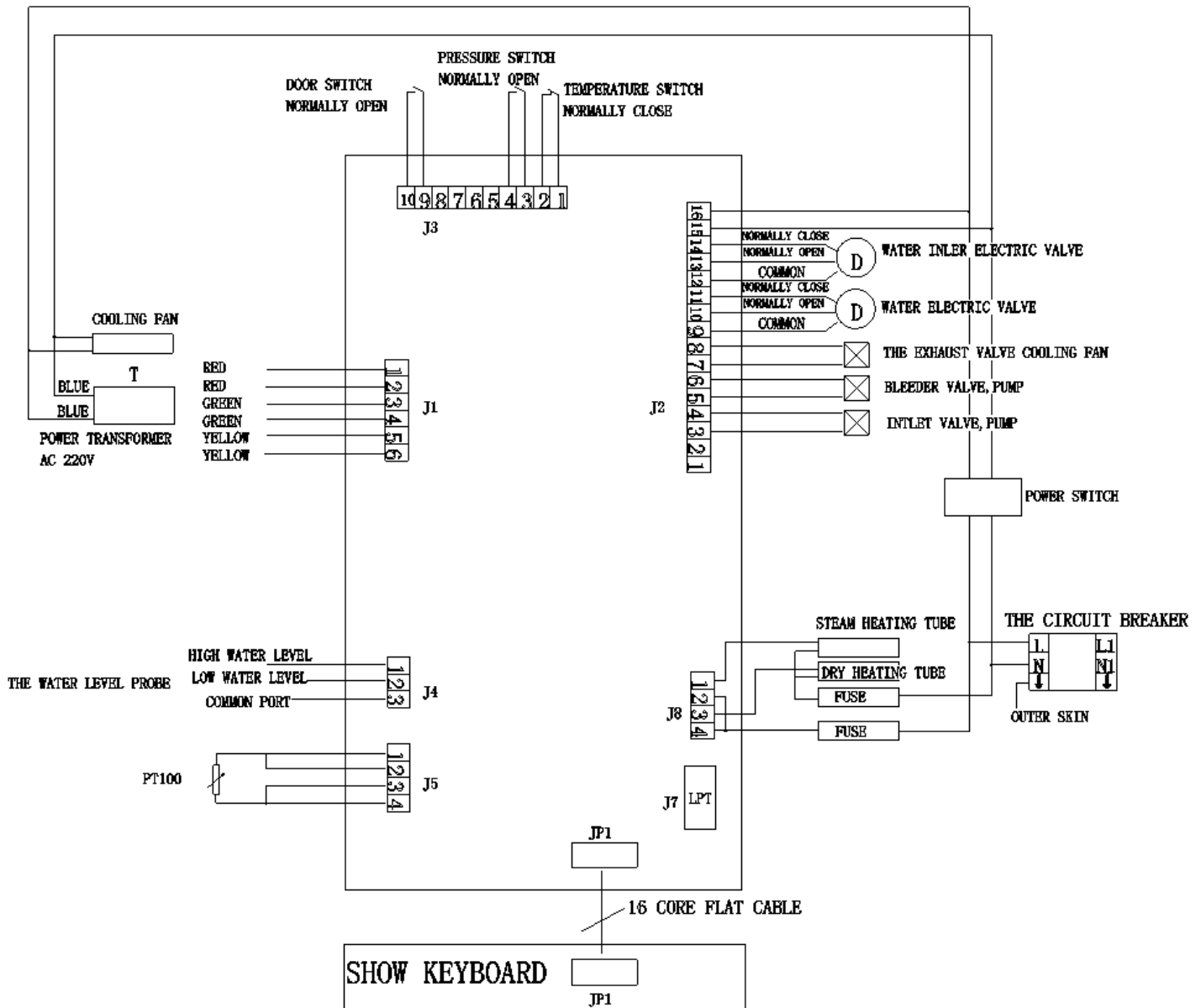

If necessary, we can provide circuit diagram and effecting the necessary information. If there is any doubt about electrical circuit repair, you can contact with our company.

06. Common Fault Analysis and Troubleshooting

| Seq | Fault Symptoms | Cause Analysis | Troubleshooting Methods |
|-----|---|--|---|
| 1 | Water level light on, temperature not rising on display | a.The sterillization time is not set b.Circuit board damaged c.Heat tube damaged | Set the sterillization time Be replaced by professional Be replaced by professional |
| 2 | Operating status, shell electrified | a.Heat tube damaged b.Power supply without grounding line. | Be replaced by professional Grounding wires by professionals. |
| 3 | Temperature display shows - - - - | a.The temperature sensor damaged b.Circuit board damaged | Be replaced by professional Be replaced by professional |
| 4 | Digital display shows - O O O or garbled characters | Circuit board damaged | Be replaced by professional |
| 5 | Pressure gauge with water vapor inside | The pressure gauge spring tube leak | Be replaced by professional |
| 6 | Pressing the shift key, no blinking in the corresponding position | The control panel damaged | Be replaced by professional |
| 7 | Display ERR—1 or door control light not on | a.The cover is not closed in place b.Light chain broken | a.Open the cover,and close the cover in place. b.Change the light chain |
| 8 | Display ERR-2 | Water level below low water level | Add enough water |
| 9 | Display ERR 3 | Sterilization temperature too high | SCR failure and the temperature sensor failure |
| 10 | Display ERR 5 | Heating element overheating protection | Temperature control switch failure (Professional repair) |
| 11 | Display ERR 6 | Pressure too high | Pressure switch, safety valve failure (Professional repair) |
| 12 | Display ERR 7 | Power outage during sterilization process | Manually unlock by pressing the function key |
| 13 | Display ERR 8 | Manual reset during sterilization process | Manually unlock by pressing the function key |



| | | | |
|----|----------------|---|--|
| 14 | Display ERR 9 | Indoor temperature sensor malfunction | Replace the sensor (to be done by a professional). |
| 15 | Display ERR 10 | Wall temperature sensor malfunction. | Replace the sensor (to be done by a professional). |
| 16 | Display ERR 12 | Heating film malfunction(for drying function). | Replace the heating film (to be done by a professional). |
| 17 | Display ERR—20 | Timeout for heating | Replace heating element |
| 18 | Display ERR—21 | Sterilization temperature lower than set temperature, temperature sensor damaged; heating element damaged | a. Replace heating element (Professional replacement) b. Replace temperature sensor (Professional repair) |
| 19 | Display ERR—30 | indicates Water inlet failure. | Check if the water source is sufficient, and ensure that the inlet pipe is unobstructed. The failure of the water pump or the inlet solenoid valve should be examined. |

PACKING LIST

| Seq | Item | Quantity | Note |
|-----|---------------------------|----------|------|
| 1 | Vertical Steam Sterilizer | 1 | |
| 2 | Operating Manual | 1 | |
| 3 | Gloves | 1 pair | |
| 4 | Bucket/Basket | 1 | |
| 5 | Air releasing tube | 1 | 2m |



LABSTAC

Email: contact@labstac.com
Website: www.labstac.com